

Direct analysis of trace elements in estuarine waters using triple quadrupole ICP-MS

Authors: Lionel Dutruch¹, Jörg Schäfer¹,
Mélina Abdou¹ and Daniel Kutscher²;
¹University of Bordeaux, Pessac, France;
²Thermo Fisher Scientific, Bremen, Germany

Keywords: Argon gas dilution,
contamination, estuarine waters,
interference removal, matrix effects

Goal

To show accurate quantification of toxic elements in a heavy sample matrix like estuarine waters using argon gas dilution.

Introduction

An estuary is an ecosystem, which is characterized by both marine (e.g. tides, or erosion through waves) and riverine processes (e.g. freshwater and sediment inputs). At the interface between fresh- and sea water domains, estuarine waters are often referred to as brackish waters with salinities between 1 and up to 35 (corresponding to total dissolved solids between 0.1–≤3.5%). Coastal zones including estuaries are historically populated by humans, hosting major cities and industrial activities. Estuaries are the focal points of aquatic contaminant transport from the



continent to the sea, facing contamination from industrial, agricultural and urban sources, especially over the past decades. Important examples for estuaries are e.g. Puget Sound in the northwestern US, Rio de la Plata in South America, or the Thames Estuary in the UK. In this study, the focus is on the Gironde Estuary in southwest France, a major fluvial-estuarine system in Europe. The main objective of this study is to track and control historical and ongoing multi-metal contamination of the Gironde Estuary waters. Clear evidence for contamination (mainly Cd) has been observed in seafood (oysters) from the estuary mouth and the nearby Marennes-Oléron Bay, Europe's major oyster production area¹.



However, the analysis of samples containing high salt loads imposes special challenges when using ICP-MS. Salts may crystallize during the process of nebulization or deposit on surfaces of the interface region, leading to blockage of nebulizers and cone orifices. This may lead to severe reduction of signals and cause drift problems. With respect to spectral interferences, most commonly chlorine based polyatomics, such as those interfering on vanadium, chromium and arsenic, need to be removed. Additionally, strong interferences can be observed on copper, based on the presence of sodium and magnesium in estuarine and sea waters.

In order to overcome the impact of the sample matrix, samples can be diluted using clean diluents, but especially for the analysis of elements at trace or ultra-trace levels, dilution always induces the risk of contamination or overdilution, leading to final concentrations in the measured solution being lower than instrumental detection limits (IDL). Another appealing alternative is the use of argon gas to dilute the sample before it enters the plasma. Although this also leads to a significant reduction in achievable instrument sensitivity, method detection limits (taking into account all steps in sample preparation) can be less compromised as compared to liquid dilution.

Method

Sample preparation

Estuarine water samples were collected in the high salinity range of the Gironde Estuary i.e. relatively close to the estuary mouth, ~100 km downstream from Bordeaux, France. Sampling has been performed onboard the research vessel *Thalia* (Ifremer) over a cycle of 30 hours, implying that the water masses sampled have variable salinity due to the strong ebb-flood cycle in this meso-/ macrotidal estuary. Salinities in the samples ranged from $S = 30.2$ to 31.8 . To avoid potential matrix effects caused by different salinity, all samples were adjusted to a salinity of $S=30$. The star in Figure 1 shows the sampling site in the Gironde Estuary mouth, being part of a larger sampling campaign along the entire estuarine salinity gradient (data not shown).



Figure 1. Sampling location along the Gironde Estuary

Due to the extremely low concentrations in seawater commonly observed for some of the analytes (especially Pb, but also Cd), careful control over potential sources of contamination and clean laboratory conditions are key to successful analysis. The labware was acid-cleaned (soaking 3 days in 10% HNO_3 Normapur®, VWR-BDH Chemicals), thoroughly rinsed with MilliQ® water (Merck), dried under a laminar flow hood in a clean lab (over-pressurized, filtered, air-conditioned atmosphere), then sealed in double plastic bags until use. All samples were filtered onboard immediately after sampling using $0.2 \mu\text{m}$ membrane filters (MINISART® NML, Sartorius), acidified (1/1,000, HNO_3 Suprapur®, Merck), and stored in the dark at 4°C pending analysis.

Instrument configuration

A Thermo Scientific™ iCAP™ TQ ICP-MS in combination with an SC-4DX Autosampler (Elemental Scientific, Omaha, NE) was used for analysis. The instrument was operated using Argon Gas Dilution (AGD) allowing direct analysis of estuarine waters without any prior dilution. Tuning of the system was accomplished using the autotune routines provided with the Thermo Scientific™ Qtegra™ Intelligent Scientific Data Solution™ Software. Typical operating conditions are summarized in Table 1. The selection of analytes, appropriate analysis mode (single quad vs. triple quad, choice of reactive gas) was accomplished using the Reaction Finder method development assistant. For the selected elements, this resulted in only arsenic being acquired in a triple quadrupole mode using oxygen as a reactive gas, whereas for all other analytes, the use of helium and kinetic energy discrimination was recommended. The mass selection in the first quadrupole was controlled using intelligent Mass Selection (iMS) in all cases.

Table 1. Instrument configuration

Parameter	Value
Nebulizer	MicroMist Quartz nebulizer 0.4 mL·min ⁻¹ , pumped at 40 rpm
Spraychamber	Quartz cyclonic spraychamber cooled at 2.7 °C
Injector	2.5 mm i.d., Quartz
Interface	High Matrix (3.5 mm) insert, Ni cones
RF power	1,550 W
Nebulizer gas flow	0.73 L·min ⁻¹
Additional gas flow (AGD)	97%
QCell settings	SQ-KED TQ-O₂
Gas flow	100% He, 4.2 mL·min ⁻¹ 100% O ₂ , 0.3 mL·min ⁻¹
CR bias	-21 V - 6.4 V
Q3 bias	-18 V -12 V
Scan settings	0.1 s dwell time per analyte, 10 sweeps, 3 main runs

General analytical condition

For calibration and quality control, a certified reference material was used (CASS 6, Nearshore Seawater Certified Reference Material for Trace Metals and other Constituents, National Research Council Canada). Matrix matched calibration curves were generated by addition of increasing concentrations of the elements investigated in this study directly into aliquots of the CASS-6 CRM. Table 2 gives an overview on the added concentrations for the different analytes. As the calibration was matrix matched no internal standard was used in the analysis. Again, all solutions were

adjusted to a salinity of S=30. Following 8–9 unknown samples, CASS 6 was repeatedly analyzed by standard addition in order to (i) check the accuracy of the method and (ii) monitor potential sensitivity drift.

Table 2. Added concentrations for calibration curves per element

	Cu	Zn	As	Cd	Pb
Zero STD	0	0	0	0	0
Standard 1 [µg·kg ⁻¹]	0.5	1.3	1.0	0.02	0.01
Standard 2 [µg·kg ⁻¹]	1.0	2.6	2.0	0.04	0.02
Standard 3 [µg·kg ⁻¹]	1.5	3.9	3.0	0.06	0.03
Standard 4 [µg·kg ⁻¹]	2.0	5.2	4.0	0.08	0.04
Standard 5 [µg·kg ⁻¹]	2.5	6.5	5.0	0.10	0.05

Results

The results obtained are summarized in Table 3. As can be seen, quantitative recoveries are obtained for all elements under study in the CASS-6 CRM. The CRM was analyzed 4 times throughout the analysis and demonstrated low relative standard deviations despite the extremely low concentrations of some of the elements. For arsenic (the only element measured in both single and triple quadrupole modes), there is no significant difference in the results considering the uncertainty information in the certificate of the CASS 6 reference material. In this sample matrix, the predominant interference on ⁷⁵As are chlorine -and calcium-based species, such as ⁴⁰Ar³⁵Cl⁺, ⁴⁰Ca³⁵Cl⁺ or ⁴⁰Ca(OH)₂H⁺. Due to their polyatomic nature, these interferences can be efficiently removed by KED alone. However, it is worth noticing that the triple quadrupole based mode using oxygen offered a much higher detection sensitivity (more than double in comparison to KED) and significantly lower detection limits (more than 5 times lower). Other interferences, such as doubly charged ions of the Rare Earth Elements, might affect the results for elements such as arsenic or selenium, and can only be removed using triple quadrupole technology. However, in this study they were not found to be causing any bias to the results.

The results of this study were compared to an earlier study using a different analytical technique, i.e. a submersible voltammetry system validated for measuring estuarine samples². The comparison shows that results are very similar for arsenic, cadmium and lead. Results for copper and zinc deviated slightly from earlier results probably due to their higher affinity for forming complexes with organic molecules present in sea water, which are not detected by in-situ voltammetry.

Table 3. Results obtained for the measurement of CASS 6 CRM and 18 samples

Element	Cu	Zn	As		Cd	Pb
Mode	SQ-KED	SQ-KED	SQ-KED	TQ-O2	SQ-KED	SQ-KED
Result CASS 6 CRM [$\mu\text{g}\cdot\text{kg}^{-1}$] (N=4)	0.57 ± 0.012	1.89 ± 0.23	1.04 ± 0.11	1.09 ± 0.08	0.027 ± 0.004	0.013 ± 0.002
Certified value [$\mu\text{g}\cdot\text{kg}^{-1}$]	0.530 ± 0.032	1.27 ± 0.18	1.04 ± 0.10		0.0217 ± 0.0018	0.0106 ± 0.0040
Concentration range in samples [$\mu\text{g}\cdot\text{kg}^{-1}$]	0.31-0.56	0.41-2.34	1.32-1.88		0.017-0.058	0.023-0.042

Conclusion

The direct analysis of estuarine waters without prior dilution is possible using AGD on the iCAP TQ ICP-MS. The results obtained for the CASS-6 CRM indicate accurate and precise quantification is possible at very low concentration levels. The results obtained for the samples collected in the high salinity range of the Gironde Estuary show that historical metal contamination in the estuarine waters persists, although at lower levels (especially for elements such as Cd or Zn) as compared to earlier studies conducted in the 1990's³. This observation fits with the continuous decrease of Cd concentrations determined in wild oysters from the Gironde Estuary mouth⁴. Comparing the data to the results of an earlier study, it is clear that both, ICP-MS and voltammetry methods, provide similar results at trace and ultra-trace levels. Slight differences in results may occur due to metal species/complexes not detected in voltammetry. If submersible voltammetry systems allow for in-situ measurements of a number of trace metals, ICP-MS is a time-efficient alternative and allows the analysis of a wider range of elements in one aspiration of the sample due to its inherent multi-elemental capability. Future work on multi-element analysis in seawater should include an even wider range of elements, including emerging metal contaminants.

References

- Dabrin, A., Schäfer, J., Blanc, G., Strady, E., Masson, M., Bossy, C., Castelle, S., Girardot, N., and Coynel, A. 2009. Improving estuarine net flux estimates for dissolved cadmium export at the annual timescale: Application to the Gironde Estuary. *Estuarine and Coastal Shelf Science* 84, 429–439.
- www.schema-ocean.eu
- Jouanneau, J.M., Boutier, B., Latouche, C., Phillips, I., 1990. Cadmium in the Gironde fluvio-estuarine system: behaviour and flow. *Science of the Total Environment* 97/98, 465–469.
- Lanceleur L., Schäfer J., Chiffolleau J.F., Audry S., Auger D., Renault S., Baudrimont M., Blanc G. (2011). Long-term (30 years) records and relationships of cadmium and silver contamination in sediment and oysters from the Gironde fluvial-estuarine continuum. *Chemosphere* 85:1299-1305.

Acknowledgement

The authors greatly acknowledge the use of the R/V Thalia (Ifremer) of the French National Coastal Research Fleet. Sampling and sample treatment was funded within the SCHeMA project (Integrated In Situ Chemical Mapping Probes; EU FP7 Ocean 2013.2; Project-Grant Agreement 614002).

Find out more at thermofisher.com/TQ-ICP-MS

For Research Use Only. Not for use in diagnostic procedures. © 2020 Thermo Fisher Scientific Inc. All rights reserved.

Normapure is a registered trademark of VWR BDH Chemicals; MilliQ and Suprapur are registered trademarks of Merck Group; Minisart is a registered trademark of Sartorius AG. SC and Elemental Scientific are trademarks of Elemental Scientific Inc. MicroMist is a trademark of Glass Expansion Ltd. All other trademarks are the property of Thermo Fisher Scientific and its subsidiaries. This information is presented as an example of the capabilities of Thermo Fisher Scientific products. It is not intended to encourage use of these products in any manners that might infringe the intellectual property rights of others. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representatives for details. Not all products are available in all countries. Please consult your local sales representatives for details. **AN44417-EN 0620C**

ThermoFisher
SCIENTIFIC