


Thermo Scientific Theta Probe

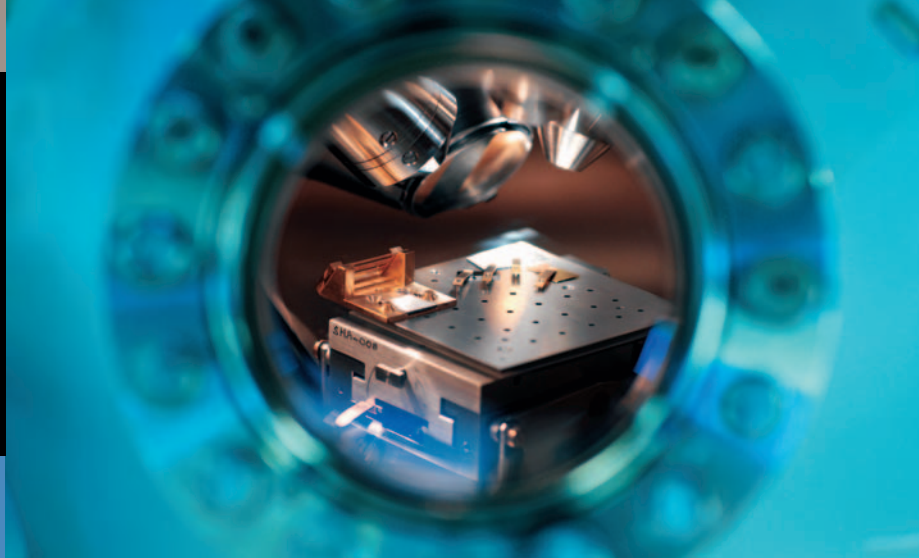

Small Spot XPS Spectrometer with Parallel ARXPS Capability



Polymers



Semiconductor



Self Assembled Monolayers



Surface Modification

Theta Probe

The Thermo Scientific Theta Probe is a major advance in surface and ultra-thin film analysis, featuring small spot XPS and a unique parallel angle-resolved XPS capability suitable for applications such as:

- Ultra-thin films
- High-k dielectrics
- Surface modification of insulators
- Self assembled monolayers
- Modified polymers

Theta Probe combines ease-of-use with rapid sample throughput.

- Fast snapshot spectra
- Maximum sensitivity in small area XPS
- Unique parallel angle resolved XPS (PARXPS) for fast data acquisition
- Chemical state and depth information
- Non-destructive depth analysis
- Overhead sample viewing for fast and accurate sample alignment
- Automated multi-sample analysis



A New Dimension in Surface Analysis

Theta Probe

Theta Probe is an exceptionally easy-to-use problem-solving tool. It delivers:

- Quantitative surface and thin film analysis
- Small area analysis with high sensitivity
- Non-destructive analysis of ultra-thin films (<10 nm)
- Sputter depth profiles
- Maps of
 - Elemental distribution
 - Chemical state distribution
 - Film thickness
 - Film uniformity

Small Area XPS (SAXPS)

Theta Probe uses a microfocusing monochromator for source-defined SAXPS, making it as simple as large area XPS

- User-selectable spot size from 15 μm to 400 μm
- Source-defined SAXPS provides maximum sensitivity especially at the smallest areas
- Moveable anode for maximum lifetime
- Effective charge compensation

Parallel Angle-resolved XPS

For advanced thin film analysis Parallel Angle Resolved XPS (PARXPS) has many advantages

- PARXPS can be used for small areas, large samples and within images and depth profiles
- PARXPS is faster than the traditional method involving sample tilting
- The analysis area remains constant, independent of electron emission angle
- The energy resolution is constant over all angles because the whole of the X-ray spot is within the field of view of the spectrometer
- Mapping can be combined with ARXPS to determine the uniformity of ultra-thin layers

Measurement

PARXPS is a powerful tool for non-destructive determination of:

- The sequence of layers in an ultra-thin film
- Thickness of the layers, including buried layers
- The concentration depth profile through the layers
- Uniformity of thickness, composition and distribution of material within the layers

Avantage





The Thermo Scientific *Avantage* data acquisition and processing software incorporates all the features necessary to extract the maximum amount of information from the analysis and has a comprehensive set of data and image processing tools including:

- Linear, Shirley and Tougaard background subtraction
- Peak identification
- Quantification
- Peak fitting
- Non-linear least squares fitting
- Multi-dimensional data display
- PARXPS tools
 - Relative depth plotting
 - Multiple overlayer thickness calculator
 - Depth profile generator

Sample Handling

Theta Probe is equipped with versatile and precise sample handling features.

- Stage movement 70 x 70 x 25 mm for large samples or a large number of small samples
- Multi-sample unattended operation for maximum productivity
- CCD camera and zoom microscope aligned with the analysis position for fast and accurate sample alignment
- Overhead optical view allows the stage position to be controlled from the optical image
- Optional 8-sample rotation holder

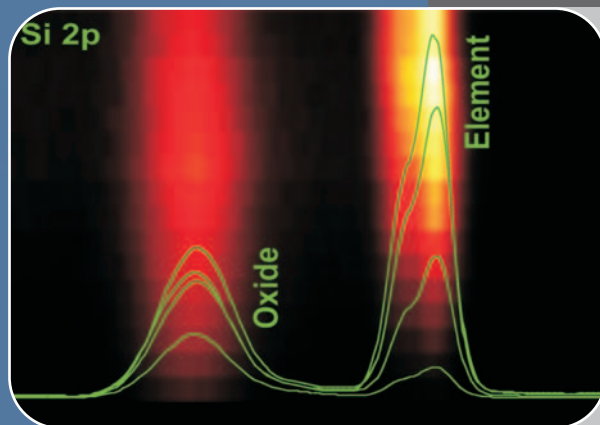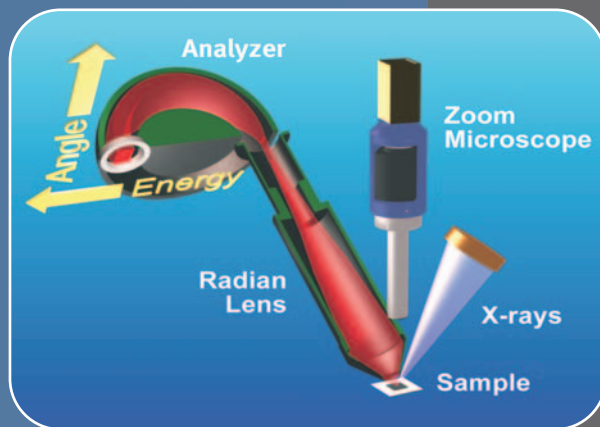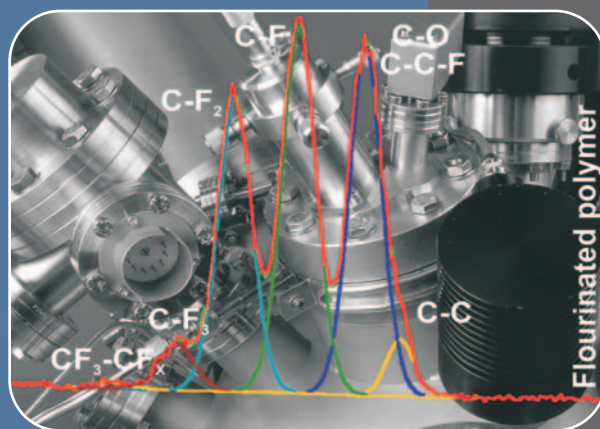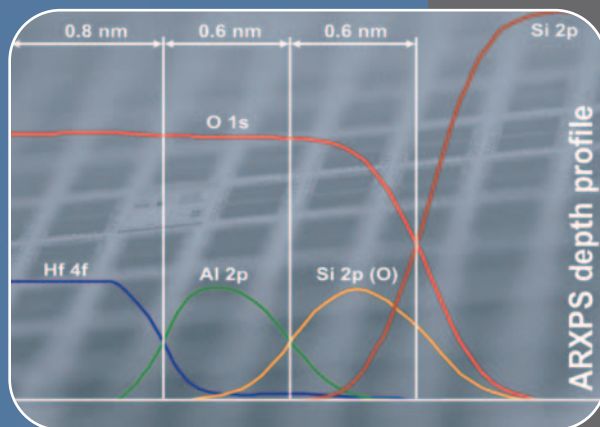

Theta Probe

Electron Analyzer

- Double-focusing full 180° spherical sector Analyzer
- Multi-element electrostatic RADIAN input lens for spectroscopy
- Two-dimensional, multi-channel spectroscopic detector for spectroscopy and PARXPS

Microfocused Monochromated X-ray Source

- 250 mm Rowland circle monochromator
- Microfocused electron gun and multi-position aluminum anode X-ray source
- Thermostatically controlled quartz crystal and crystal manipulator
- Spot sizes down to 15 µm

Combined Low-energy Electron/Ion Flood Gun

- Charge neutralization
- Flood gun assembly incorporating electrostatic deflection for precise alignment
- Low energy-spread, high brightness electron source

Ion Gun

- Sample cleaning
- Depth profiling
- Beam scanning and video
- Differential pumping

Sample Viewing


- X-ray shielded viewport
- Zoom microscope sample alignment facility with CCD camera
- Multiple light sources

Fast Entry Chamber

- UHV compatible sample transfer mechanism
- Fully safety interlocked pneumatic 98 mm gate valve to analysis chamber
- Turbomolecular pump and rotary backing pump

5-axis Sample Stage

- High precision, automated sample stage with internal stepper motors
- Stage controller with trackball and joystick control interfaced to the *Avantage* data system
- Maximum analysis area 70 mm x 70 mm
- Maximum sample thickness 25 mm

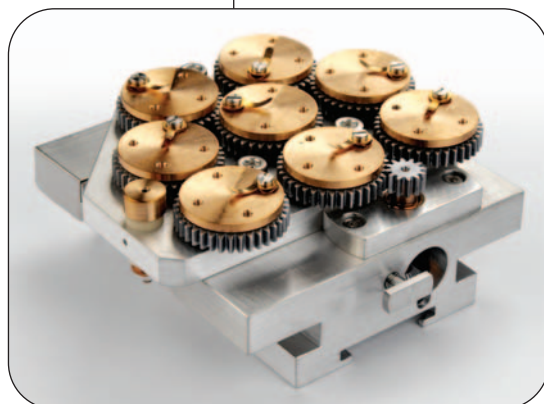

In addition to these offices, Thermo Fisher Scientific maintains a network of representative organizations throughout the world.

Africa

+43 1 333 5034 127 • analyze.emea@thermo.com

Australia

+61 2 8844 9500 • analyze.au@thermo.com

Austria

+43 1 333 50340 • analyze.at@thermo.com

Belgium

+32 2 482 30 30 • analyze.be@thermo.com

Canada

+1 800 530 8447 • analyze.ca@thermo.com

China

+86 10 8419 3588 • analyze.cn@thermo.com

Denmark

+45 70 23 62 60 • analyze.dk@thermo.com

Europe - Other

+43 1 333 5034 127 • analyze.emea@thermo.com

France

+33 1 60 92 48 00 • analyze.fr@thermo.com

Germany

+49 6103 408 1014 • analyze.de@thermo.com

India

+91 22 6742 9434 • analyze.in@thermo.com

Italy

+39 02 950 591 • analyze.it@thermo.com

Japan

+81 45 453 9100 • analyze.jp@thermo.com

Latin America

+1 608 276 5659 • analyze.la@thermo.com

Middle East

+43 1 333 5034 127 • analyze.emea@thermo.com

Netherlands

+31 76 579 55 55 • analyze.nl@thermo.com

South Africa

+27 11 570 1840 • analyze.sa@thermo.com

Spain

+34 914 845 965 • analyze.es@thermo.com

Sweden / Norway / Finland

+46 8 556 468 00 • analyze.se@thermo.com

Switzerland

+41 61 48784 00 • analyze.ch@thermo.com

UK

+44 1442 233555 • analyze.uk@thermo.com

USA

+1 800 532 4752 • analyze.us@thermo.com

www.thermo.com



VG Systems Ltd. Trading as Thermo Fisher Scientific, East Grinstead, UK is ISO Certified.

©2008 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.

Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

BR31003_E 05/08M